

Lindberg Blue M Convection Oven

Convection Oven	
	
Tool Type: Oven	
Location: Microfluidics Lab	
Description: Mechanical Convection Oven	
Manufacturer: Lindberg / Blue M	

About

Safety Concerns

High Voltage, High Temperature, no Combustibles or Hazardous Materials

Operating Procedures

1. Enter the set point temperature (see instructions for the different controllers below)
2. Place your items in the oven. Remember to place liquid materials on a baking sheet or other secondary container.
3. Set the countdown timer for the length of your run. If the timer reads zero time remaining another user who needs to use the oven may remove your items.
4. Remove your items from the oven and close the door.
5. The oven is left powered on.

Detailed Specifications

Reference Documentation

[Laboratory Oven SOP](#)

Operations Manual

From: <https://microfluidics.cnsi.ucsb.edu/wiki/> - **Innovation Workshop Wiki**

Permanent link: <https://microfluidics.cnsi.ucsb.edu/wiki/doku.php?id=lindberg-convection-oven&rev=1672952693>

Last update: **2023/01/05 21:04**

